

Underivatized Amino Acids

Overview

An introduction is needed as to which unit these results are from.

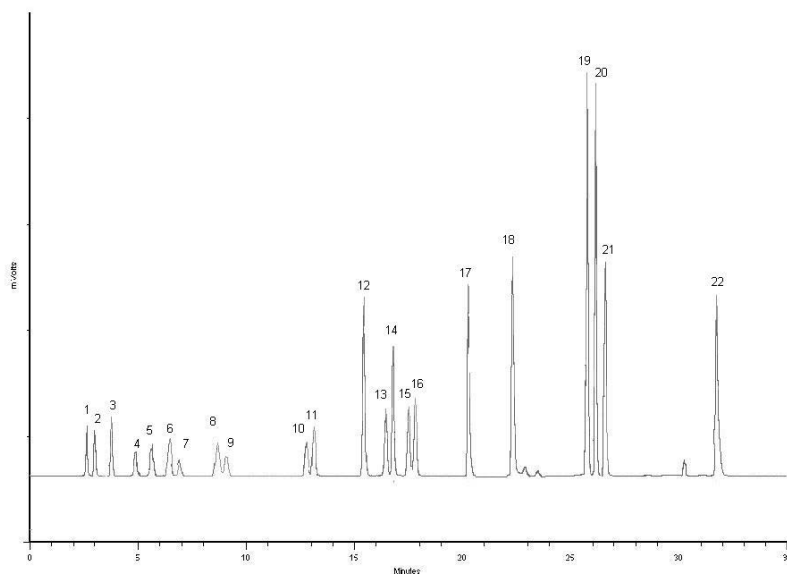


Figure 1: HyperCarb Column

- Mobile Phase: A: Nonafluoropentanoic acid B: ACN/ Nonafluoropentanoic acid
- Sample: 10nM Each Amino Acid
- Peaks: 1:Gly, 2: Ser 3: Ala, 4: Thr, 5: Asn, 6: Pro, 7: Asp 8: Gaba, 9: Gln, 10: E, 11: V, 12: Lys, 13: Leu, 14: Cys, 15: Met, 16: Ile, 17: His, 18: Arg, 19: Phe, 20: Tyr, 21: Impurity, 22: Trp
- Spray Chamber: 30°C
- Drift Tube: 55°C
- Filter: 6

Teledyne Isco

P.O. Box 82531, Lincoln, Nebraska, 68501 USA
Toll-free: (800) 228-4373 • Phone: (402) 464-0231 • Fax: (402) 465-3091
E-mail: IscoInfo@teledyne.com

Teledyne Isco is continually improving its products and reserves the right to change product specifications, replacement parts, schematics, and instructions without notice.