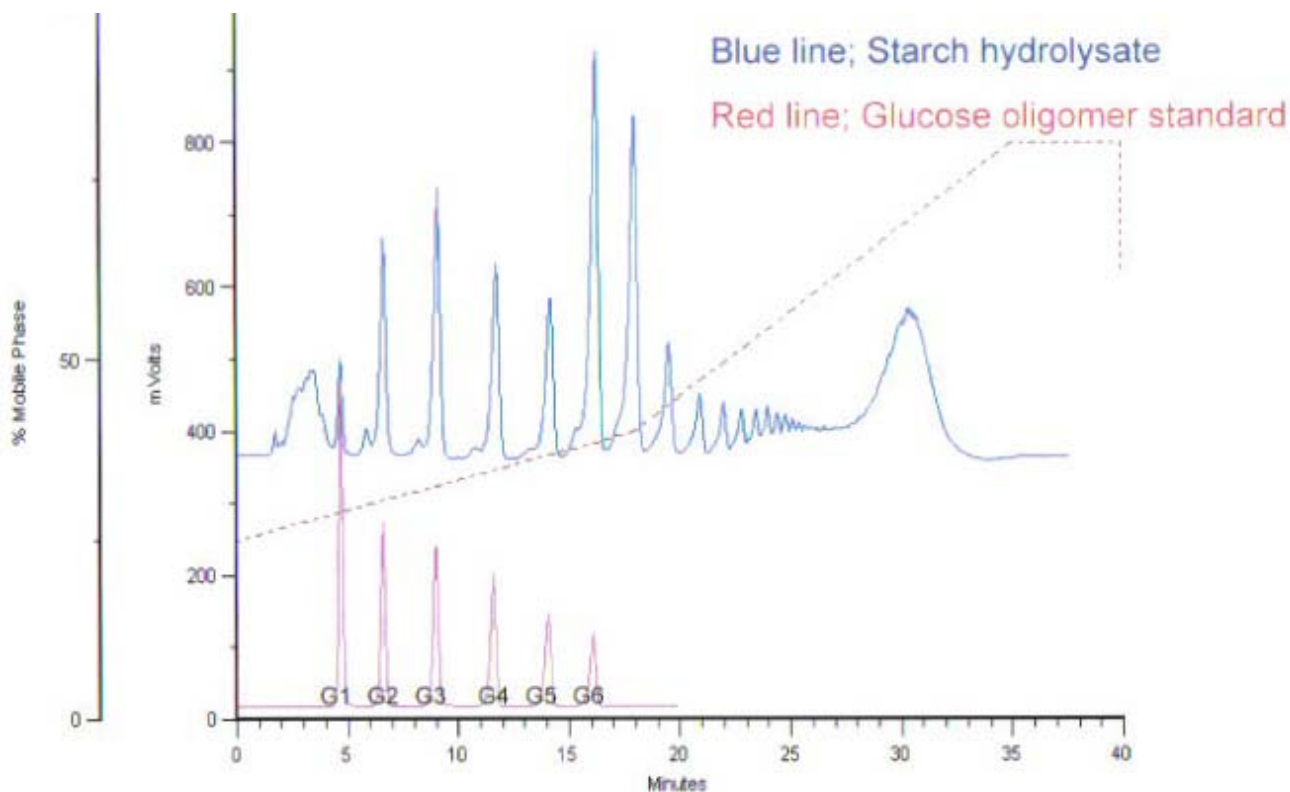


# Starch Hydrolysate and Glucose Oligomer

## Overview

Need an overview



**Figure 1: need a caption**

- Column: Asahipak NH2P 4.6mm x 150mm
- Column Temperature: 40 °C
- Mobile Phase: A: CH<sub>3</sub>CN B: Water
- Gradient: 0 min 25% B, 18 min 40% B, 35 Min 80% B
- Flowrate: 1ml/min
- ELSD: Model 200
- Spray Chamber Temperature: 25 °C
- Drift Tube Temperature: 45 °C
- Gain: Normal
- Full Scale: 5V

## Teledyne Isco

P.O. Box 82531, Lincoln, Nebraska, 68501 USA

Toll-free: (800) 228-4373 • Phone: (402) 464-0231 • Fax: (402) 465-3091

E-mail: [IscoInfo@teledyne.com](mailto:IscoInfo@teledyne.com)

Teledyne Isco is continually improving its products and reserves the right to change product specifications, replacement parts, schematics, and instructions without notice.